## **30" Stainless Ductless Fume Hood**

## Model # SS-330-DCH-ST

MADE IN THE USA

Since 1983

The **30" Wide Stainless Ductless Fume Hood** combines powerful air flow with a sturdy and compact enclosure for optimal respiratory and environmental protection. This fume hood helps protect the operator and environment from hazardous fumes or moderately corrosive particulate generated from applications performed within the hood.

The **30" Stainless Ductless Fume Hood's** powerful fan directs fumes into the filter chamber containing high quality filter media selected depending on the application. Filter media choices include **HEPA** [up to 99.97% efficiency on particles 0.3 microns and larger], **ASHRAE** [up to 95% efficiency on particles 0.5 microns and larger, **Activated Carbon**, and **specialty-blended filtration** [i.e. acid gas, mercury, aldehyde, and ammonia].

Typical uses for the **30**" **Stainless Ductless Fume Hood** include chemical fumes, pharmaceutical compounding, soldering applications, dust removal, biological applications, and a variety of applications requiring fume and dust removal. The **30**" **Stainless Ductless Fume Hood** can provide an effective and economical solution for many commercial, medical, laboratory and industrial applications that require a small containment area. For additional visibility, a transparent version is available at extra cost.

Our line of ductless fume hoods is not intended for sterile compounding applications.



30" Stainless Ductless Fume Hood



30" DCH provides fume control for a various applications.

## **Product Features**

- Includes: variable speed controller (115V only), velometer, and LED light
- Optional: clear anti-static vinyl curtains,
   Magnehelic Gauge, hour counter, and spill tray
- **Standard Sizes:** 30", 40", and 50" (custom sizes available)

- · Quiet operation
- Long filter life
- · Simple, quick "no tool" filter change
- Portable and lightweight
- Reliable, low maintenance operation



## **Product Specifications**





Shown with outer dimensions. Dimensions are approximate. Shown with optional curtains.

Hood Dimensions (D x W x H)	Outer: 24" x 32" x 25" Interior: 23.5" x 30" x 24" Access Area: 8" x 30"
Hood Material	Hood: 16 awg 304 stainles steel Fasteners: 304 stainless steel Front Window: 0.25" clear polycarbonate
<b>Blower Dimensions</b>	15.5"L x 13" W x 14.5" H
Blower Material	16 Ga. carbon steel with polyester epoxy coating on all surfaces
Weight	100 lbs
Air Volume	350 CFM
(approximately)	(varies with filter media)
Inlet Velocities on High (approximately)	Carbon Pre-filter + HEPA Filter - 130 FPM high, 25 FPM low Carbon Pre-filter + Heavy Duty Carbon Filter - 115 FPM high, 20 FPM low
Electrical	115/1/60, 2 amps 220/1/50, 0.75 amps 8' grounded power cord with NEMA 5-15P plug
Sound Level	61 to 66 dba (low to high)
Maximum Temperature	Operating temperature not to exceed 250° F
Main Filter Options	HEPA (up to 99.97% efficient on particles down to 0.3 microns)  ASHRAE (up to 95% efficient on particles down to 0.5 microns)  Activated Carbon  Specialty-blended filter media (i.e. acid gas, mercury, aldehyde, ammonia)  Limited two-year warranty
Warranty	from date of shipment on defects due to materials or workmanship.
Patent #	5,843,197



21221 FM 529 Rd Cypress, TX 77433 1.800.799.4609 sales@sentryair.com

www.sentryair.com