

## Walk-in Hood Control Panel

Model # SS-WICP


The **Walk-in Hood Control Panel** provides a central area to control the air cleaners and LED lights and a digital display that helps the operator monitor the enclosure. The control panel hardwires into the system using heavy-duty umbilical cords and mounts on the side of the structure. The **Walk-in Control Panel** includes a *master power switch* to all the fan units, *adjustable LED light controls*, and a *digital display showing the digital Magnehelic gauge and hour counter*. The *digital Magnehelic gauge* measures the static pressure of each unit's filter and helps aid as a tool to inform the operator when to change the filters. The static pressure will rise, and once it reaches a certain level, it indicates it is time for a filter change. The digital display rotates through each unit's pressure. The *digital hour counter* helps keep consistent maintenance schedules and can be reset after a filter change. This control panel can only be used with Model 300 and 400 fan units with up to 4 units. It does not retrofit to currently installed walk-in enclosures.

The **Walk-in Hood Control Panel** offers several safety features to help protect the operator. The control box can only be serviced by a licensed electrician and the system will turn OFF if the control panel door is opened. Also, the **Walk-in Hood Control Panel** is UL listed, under UL standard 61010-1, providing vital electrical safety.

***This easy-to-use control panel helps simplify using and operating a Sentry Air Walk-in Hood.***

### Product Features


- Connects with 1-4 Model 300 or Model 400 Air Cleaners
- Master power switch for fans
- Adjustable LED light controls. 1 LED light per air cleaner. Provides up to 2884 lumens per light.
- Digital display with digital magnehelic gauge (*for each unit*) and hour counter
- For use with new walk-in enclosure units only (*does not retrofit*)
- Safety mechanism turns OFF system when control panel door is opened
- UL Listed - Standard 61010-1 , File #: E466130
- Made in the USA

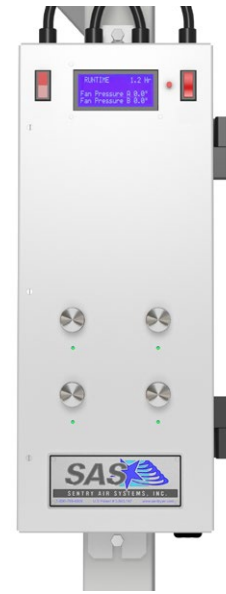

Walk-in Hood Control Panel



Walk-in Control Panel Digital Display

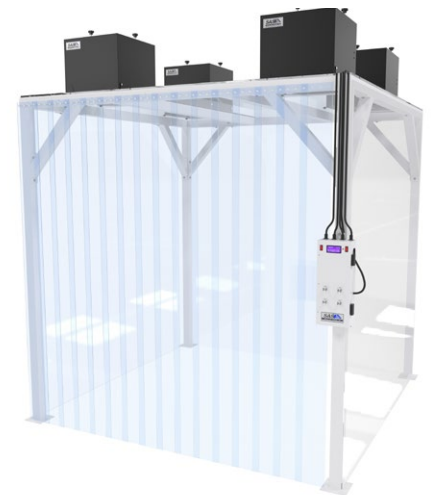

Walk-in Control Panel mounts on the side of the Walk-in Hood.




21221 FM 529 Rd, Cypress, TX 77433  
1.800.799.4609 • sales@sentryair.com

[www.sentryair.com](http://www.sentryair.com)

# Walk-in Fume Hood

## Product Specifications

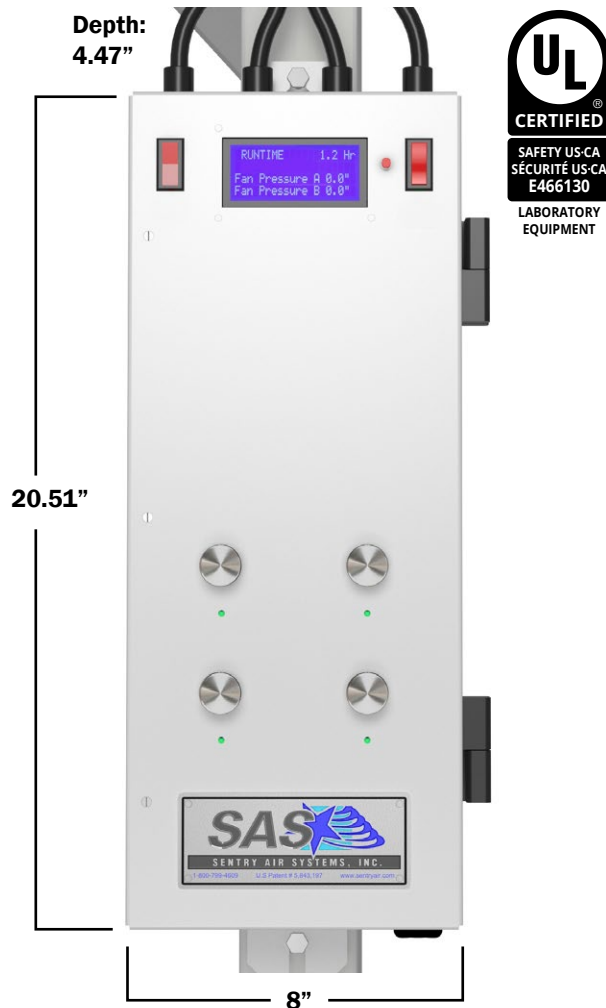

Shown with outer dimensions.  
Dimensions are approximate.



21221 FM 529 Rd  
Cypress, TX 77433  
1.800.799.4609  
sales@sentryair.com

[www.sentryair.com](http://www.sentryair.com)

<b>Panel Dimensions (D x W x H)</b>	4.47" x 8" x 20.51" (22.75" with mounting hardware)	
<b>Cabinet Material</b>	0.093" powder-coated aircraft aluminum	
<b>Weight (approximately)</b>	10.5 lbs	
<b>Electrical (Amps shown for Model 400 Units)</b>	<b>1 unit</b>	115/1/60, 2.7 amps
	<b>2 units</b>	115/1/60, 5.3 amps
	<b>3 units</b>	115/1/60, 7.9 amps
	<b>4 units</b>	115/1/60, 10.5 amps
<b>Max Wattage</b>	1188 watts	
<b>LED Light</b>	1 LED light per air cleaner 26.28" L CWH12x60-24V 2884 lumens per light	
<b>UL Listed</b>	<b>Standard:</b> 61010-1 <b>File #:</b> E466130	
<b>Disconnect Methods</b>	External Branch Circuit Breaker	
<b>Tightening Torque</b>	12 in lbs.	
<b>Warranty</b>	Limited one-year warranty from date of shipment on defects due to materials or workmanship.	